CL6-H Enclosed Standard of Spectral Irradiance User notes




Overview

The Bentham CL6-H is an enclosed spectral irradiance standard. calibrated at a distance of 5.5mm with respect to the front face of the device. This calibration is traceable to the NPL. Please refer to the accompanying certificate for calibration results.

Interchangeable adaptors to the Bentham range of input optics are available to ensure location at the calibration distance.

Mechanical

The CL6-H should be operated in the horizontal position. The appropriate DAR diffuser adaptor ring can be screwed into the front face plate, to which can be connected the input optic for system calibration.

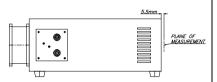

After lamp operation, allow the lamp to cool down for a few minutes prior to moving, to prevent damage to the filament, and never run the lamp with the cover in place.

Lamp Operation

Calibration distance between optic/ CL6

It is recommended to operate the lamp from a constant current supply such as the Bentham 605/608. The following provide a good guide to correct lamp operation, and to preserve the validity of the calibration.

•Ensure the correct polarity is respected at all times

- •This lamp has been calibrated with the pins down, lamp vertical, and should be used in this way. Operation in any other sense may invalidate calibration and curtail lamp life
- •Do not remove bulb from holder

•Do not touch bulb with bare fingers

•Bentham recommend that the device should be re-calibrated every 100hrs use or 1 year. whichever comes first

- •Do not run the lamp at a current lower than that at which it was calibrated
- •The lamp requires approximately five minutes warm-up time

Furthermore, it is of use to note the voltage displayed on the 605/ 608 LCD. Whilst this is for indication only, it can be used to determine lamp condition. These lamps tend to fail via either catastrophic rupture or bridging of the filament. In the latter case, adjacent filament struts collapse, short-circuiting part of the filament: at a given current the lamp voltage is therefore reduced.

Source	Current Supply (A)	Typical Operating Voltage (V)
150W	6.300	21.5

Bentham Instruments Limited 2 Boulton Road, Reading, Berkshire, RG2 0NH, UK Tel: +44 (0)118 975 1355 Fax: +44 (0)118 931 2971 Email: sales@bentham.co.uk Internet: www.bentham.co.uk

The Quartz Halogen Lamp


During use, an incandescent lamp filament reaches very high temperatures, around 3000K, at which temperature some of the tungsten, of which the filament is composed, evaporates and moves around inside the bulb by convection. Should the envelope be cold, it is likely that the tungsten is deposited thereupon, leading to gradual discolouration, as often seen in such household lamps.

Furthermore, as the tungsten is evaporated, the structure is of the filament is gradually thinned in places, leading to runaway failure: as the resistivity of the thinned sections increase, so does power dissipation and temperature, with as consequence a higher rate of evaporation. These factors conspire to thin the filament further until such point that the filament breaks, or "bridges".

The guartz halogen lamp was introduced to mitigate these problems with two main design changes brought to promote what is termed the halogen cycle.

Firstly, in order to maintain the envelope at high temperature, it is brought as close to the filament as possible (without melting). Quartz, which can withstand higher temperatures than conventional glass, is used as the envelope material.

Furthermore, the bulb is filled with a halogen gas which combines with the evaporated tungsten from the filament. Provided that the envelope of the bulb is sufficiently hot, the tungsten-halogen compound does not condense thereupon, and eventually returns via convection to the filament, at which

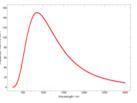

Typical spectral irradiance of CL6-H

point the tungsten is re-incorporated, releasing the halogen to continue the process.

Envelope temperature being important for the halogen cycle to operate, the lamp must be run for long enough to heat up the lamp, and should not be under-run. Operating the lamp at less than 80% of rated voltage will result in failure of the halogen cycle due to the lower bulb temperature.

Furthermore, the orientation of the lamp may become important where the envelope is not equidistant above/ below the filament. Running such a lamp in the pins down orientation is satisfactory since the upper part remains hot through convection, but in the pins down orientation, it is possible that it becomes sufficiently cool to condense tungsten. It is possible to re-evaporate tungsten deposited on the envelope by letting the lamp run at a level at which shall re-initiate the halogen cycle.

The lamp operating conditions provided overleaf lead to moderate under-running along with which comes an increase in lamp lifetime to a few thousand hours.

Finally, whilst the emission of the tungsten filament is close to that of a black body, it should be noted that it is modified by the transmission of the envelope material (without which the filament would quickly oxidise and fail). Quartz transmits over the region 200-3000nm with an absorption feature around 2500nm.

WEEE statement:

Bentham are fully WEEE compliant, registration number is WEE/CB0003ZR. Should you need to dispose of our equipment please telephone 0113 385 4352 or 4356 quoting account number 135419.



Bentham Instruments Limited 2 Boulton Road, Reading, Berkshire, RG2 0NH, UK Tel: +44 (0)118 975 1355 Fax: +44 (0)118 931 2971 Email: sales@bentham.co.uk Internet: www.bentham.co.uk

