608 Programmable Constant Current Power Supply User Notes



Overview

The Bentham 608 is a programmable constant current power supply designed for use with stabilised light sources and calibration standard sources.

Max output current= 12A dc. max output power = 250W, max compliance voltage= 35V dc

Current may be set either via a front panel keypad or via the USB interface. A socket for lamp fan operation is situated on the lower right of the front panel.

Specifications

30 ppm/℃
Green LED= Power on
Red LED = Lamp failure
Output voltage indicator
(accuracy ±0.1V- not precision meter)
110V ac (typical 200mA)
Back panel switch-
select 110V or 220/240-CV AC
50 or 60 Hz
220/240V -
2.5A anti surge
110V -5A anti surge

Lamp Operation

- Type in the required current
- The operating conditions of the various lamps supplied by Bentham are detailed overleaf
- Should an erroneous value be input hit cancel
- Hit enter/ confirm to power on lamp
- To prolong lamp life, ensure lamp polarity respected, and that those housings requiring a fan have the fan connected into the 605 lower right socket.
- Hit output off to extinguish lamp

Calibration Sources:-

· .			
Source	Current (A)	Nominal Voltage (V)	Fan
CL2	8.500	12	N
CL2-250	10.400	24	N
CL2-Nernst	8.500	4	N
CL6	6.300	24	Υ
SRS8	4.000	12	N
SRS12	8.500	12	N
TSRF/TLF_20W	3.200	6	N
TSRF/TLF1000	8.500	12	N
TSRF/TLF1800	10.400	24	N

General Illuminators:-

Source	Current (A)	Nominal Voltage	Fan
		(V)	
IL1	8.500	12	Υ
IL10	8.500	12	Y
IL6	8.500	12	Y
IL6_250	10.400	24	Υ
IL6_Nernst	8.500	4	Υ
IL7_Xe	8.500	-	Υ
IL75 Xe	5.400	-	N

Operation Notes

For correct lamp operation (and in the case of calibrated sources, to preserve the validity of calibration), please respect the following:-

All sources:-

- •Ensure the correct polarity is respected at all times, connecting red to red and black to black from lamp to 605.
- •Do not touch bulb with bare fingers.
- •Do not run the lamp at a current lower than that at which it is specified
- •The lamp should be operated in configuration in which calibrated ie. pins up/ pins down. Operation in any other sense may invalidate calibration and curtail lamp life.
- •In general bare lamps require around a 5minute warm up time, housed lamp up to 15 minutes for optimal stability
- •It is of use to note the voltage displayed on the 605 LCD. This is for indication only, but can be used to determine lamp condition1. These lamps are operated at slightly under their nominal rating, and as such the voltage readings may be lower than nominal voltage

Calibrated sources:-

- Do not remove bulb from holder.
- •Bentham recommend that the device should be re-calibrated every 100hrs use or 1-year, whichever comes first.

USB control

Please refer to Bentham USB instruments manual for generic USB control instructions.

Setting up Mains Voltage

Whilst the mains voltage is set up at Bentham according to where the device is sent, it is of good practice to verify before powering up the 605.

A toggle selector should be found underneath the mains connector to the rear of the 605. Please ensure that the voltage of the country in which the 605 is to be used is displayed.



At Bentham, the appropriate fuse (plus a supplement) is fitted. Should the voltage setting be wrong, please ensure that the correct fuse is installed.

WEEE statement:

Bentham are fully WEEE compliant, registration number is WEE/CB0003ZR. Should you need to dispose of our equipment please telephone 0113 385 4352 or 4356 quoting account number 135419.



Bentham Instruments Limited 2 Boulton Road, Reading, Berkshire, RG2 0NH, UK Tel: +44 (0)118 975 1355 Fax: +44 (0)118 931 2971 Email: sales@bentham.co.uk Internet: www.bentham.co.uk

¹⁰ne of the failure mechanisms of such lamps is "bridging" or short-circuit of part of the filament, leading to a correspondingly lower voltage and reduced light output.